
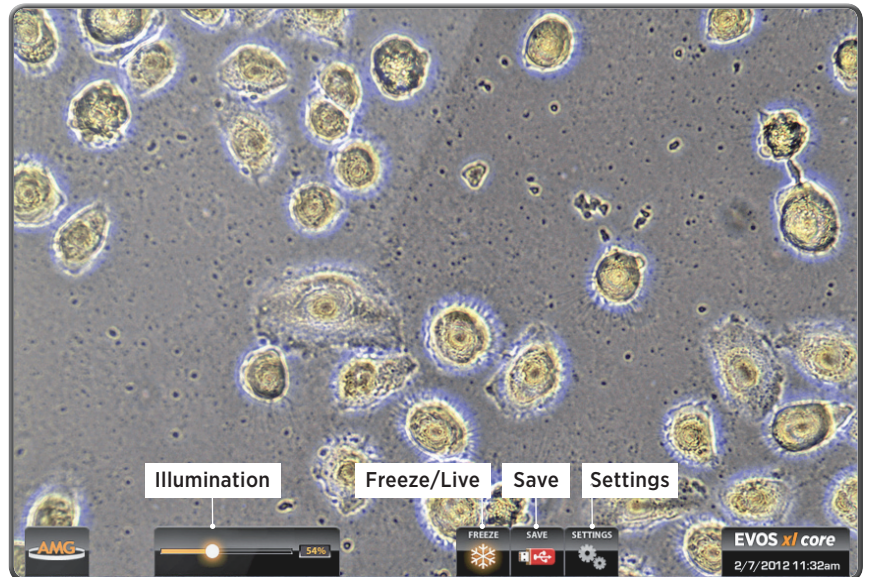


QUICK START GUIDE

- Unpack the **glass stage plate** and insert it into the **stage opening**.
- Startup takes approximately one minute. If the display remains black after startup, adjust the **illumination**.
- Set the **phase turret** position to match the selected objective and contrast method.
- The **Freeze button** toggles between freeze frame and live viewing.
- The **EVOS xl core User Guide PDF** is on the USB flash drive in the box. You can also download it from the EVOS xl core product page at www.amgmicro.com.

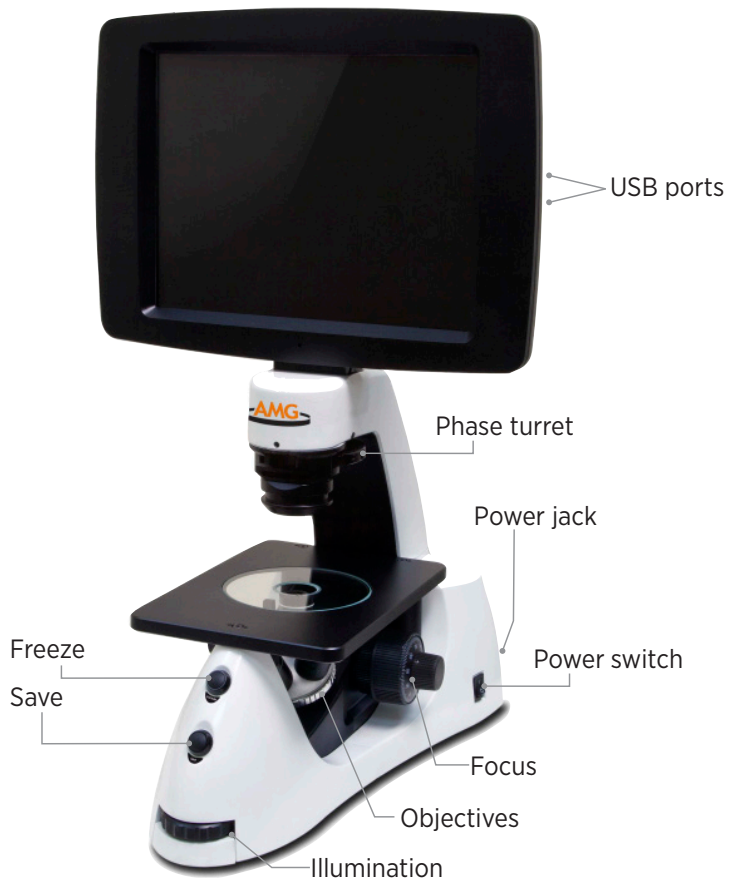

EVOS[®] xl core
Digital Inverted Microscope